

# Measurement of DHBT(Diethylhexyl Butamido triazone)

[SJ-MM-023]

2011-08-30

Version 1.0

Supersedes prior issues

Quality Assurance

## Test procedure

1. Put the sample to 10ml volume flask.
2. Add 10ml of methanol and sonicate the flask for 1hr.
3. Centrifuge (3,500rpm, 10min) & Syringe filter (0.45um) to remove impurities.

## 4. Analysis

## 5. Analysis condition

Apparatus Model	Agilent 1200 HPLC
Column	ZORBAX Eclipse XDB-C18 250 × 4.6mm I.D. 5um ZORBAX Eclipse XDB-C18 12.5 × 4.6mm I.D. 5um
Mobile phase	0.1% TFA in water : MeOH = 2 : 98
Detector	UV 300 nm
Column oven temp	35 °C
Flow rate	0.6ml/min
Injection volume	1ul



HPLC : High Performance Liquid Chromatography